

Screen 2: Automatic operation menu

Auto Operation		
Run-A	Wet gas Cycle	0
Start	Dry gas cycle	0
Reset		Return

It contains 4 buttons and 2 feedback variable status boxes, automatic operation will start when Run-A and Start are pressed (press Run-A first then Start), Wet gas Cycle box will now display wet gas clean cycle value; Dry gas Cycle box will now display dry gas clean cycle value; When two cycle is completed, 1 will be incremented in the box. When all cycles are done until the set value is reached, the counting will stop and vacuum pump will keeps running. When vacuum degree reaches the set value (P3), the system will give an alarm which indicates that canisters have been thoroughly cleaned and close the valve of the canisters. Now press “Return” to return to main menu.