

Screen3: Manual operation menu

Manual Operation		
Run-M	P1	OFF
Pump	P2	OFF
Fill wet gas	P3	OFF
Fill Dry gas		Reset
		Return

It contains 6 buttons and 3 feedback variable status boxes. Press “Run-M” to start manually operation and then continue other necessary operations. “Reset” is used to reset the operation and “Return” is to return to main menu.

Screen 4: Parameter modification menu

Parameter Set	
Vent Off-Delay	100
Wet Gas Cycle	3
Dry Gas Cycle	2
Return	

It contains 1 timing units Vent Off-Delay and 2 counting units Wet Gas Cycle, Dry Gas Cycle. The value in the display box is default value, press “Return” is to return to main menu. In case data in a certain