

2. Avoid intense vibration when the system is working. Do not put stuff on the machine and the top, so that the heat can easily scatter.
3. The main equipments should put aside from strong magnetic field. The good ground is necessary for instrument operating.
4. It is forbidden to pull out electrified socket or wires so that inner apparatus will not be damaged.
5. It is forbidden to change PLC parameters without permit. If change is needed, it only can be done by specialists.
6. Ensure the plug in power for vacuum pump is 220VAC, it is forbidden to use 110 VAC power.
7. Please refer to attached “Instructions for use of vacuum pump” for maintenance and common debugs for vacuum pump.