


3.0 Installing the batteries

This unit uses 2 AA alkaline batteries.

To install batteries, extend the wire stand. Open the cover. Insert the batteries with the polarity (⊕ and ⊖) correctly aligned. Close the cover. (Figures 2 and 3)

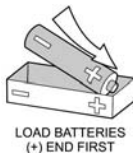

Precautions for battery replacement:

- Load the new batteries with their polarity (⊕ and ⊖) aligned correctly.
- Do not use rechargeable batteries.

4.0 Battery power consumption

4.1 Battery lifetime

The battery lifetime is dependant on the number of options the user has enabled.

The unit is shipped with the most energy demanding options disabled (Table II).

The power saving functions can be changed.

- See Section 8.2: Adjust LCD Character Contrast (p.9).
- See Section 8.3: USB Activation (p.10).
- See Section 8.4: Adjust LCD Image Backlight (p.10).
- See Section 8.7: Adjust Auto Shutoff Duration (p.12).

4.2 Battery charge indicator

The unit includes a battery charge indicator. Replace batteries as needed.

- See Section 8.5: Show Battery Charge Indicator (p.11).

Figure 2 Extend the wire stand before opening battery door.



Figure 3
Insert the batteries as marked.

Table II Default settings for the ProFLOW 6000

Auto shutoff duration	6 minutes
LCD backlight	0 (off)
LCD character contrast	5
USB port	disabled

NOTE: Store your ProFLOW 6000 in its protective storage case following use. Keep the manual under the unit; placing the manual on top can result in the unit being turned on when the lid is closed.