

If you do not already have serial terminal software, free, open source options are available online (i.e. RealTerm, etc.). You can download RealTerm\* software from:

<http://realterm.sourceforge.net/>

After installation, click the Port tab and set the following:

Baud: 57600

Port: the appropriate VCP for your ProFLOW

Parity: None

Data Bits: 8 bits

Stop Bits: 1 bit

HardwareFlow Control: None

Go to the USB menu entry in the ProFLOW and turn the transmission on. Finally, click **Open** on the Port tab in RealTerm and you will see the serial data stream begin in the terminal window (Figure 8).

If you would like to log the flow data, this can be done by clicking on the Capture tab. Set File to the name and location of the log file that you would like to save and click either the **Start: Overwrite** or **Start: Append** buttons appropriately (Figure 9).

*\*This software is not supplied or supported by Restek. User assumes all responsibility for the downloading and use of the program.*

**Figure 9** Screen capture of data collection.

