

8.0 ProFLOW 6000 menu structure


8.1 Unit power up/power down


Press the  (power) button:

The LCD screen will display the device intro screen:



Followed by measured flow data:

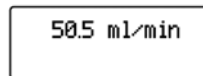

To power off, press and hold the  (power) button:

Other messages encountered at power up

Calibration
-expired-


Appears temporarily if the unit's calibration has expired (>1 year).



→ See Section 14.0: Calibration and Service


under range

If the unit is hooked up to a flow stream with a flow rate less than 0.50 mL/min. the unit will report an “under range” status for the flow. This message will appear until the flow rate exceeds 0.50 mL/min.



8.2 Adjust LCD character contrast


Press the  (menu) button.

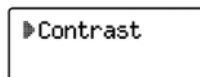
Use the   (arrow) keys to select the contrast menu:



Press the  (menu) button again to enter the value select screen.

Use the   (arrow) keys to select the contrast value.
Contrast Values: 1 (lightest) and 5 (darkest).

To return to the main menu screen, press the  (menu) button.

To exit and return to measuring flow, press the  (menu) button again.