


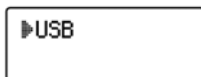




8.3 USB activation


To enable the USB, press the  (menu) button.

Use the   (arrow) keys to select the USB menu:



Press the  (menu) button again to enter the value select screen.


Use the   (arrow) keys to toggle between USB 'on' and 'off'.



To return to the main menu screen, press the  (menu) button.

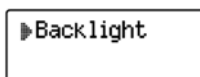
To exit and return to measuring flow, press the  (menu) button again.

→ See Section 7.0: Data Collection on the PC



8.4 Adjust LCD image backlight


Press the  (menu) button.


Use the   (arrow) keys to select the backlight menu:



Press the  (menu) button again to enter the value select screen.

Use the   (arrow) keys to select the backlight value.
Backlight Values: 0 (off) and 5 (maximum).

To return to the main menu screen, press the  (menu) button.

To exit and return to measuring flow, press the  (menu) button again.