


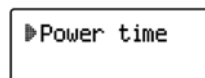




## 8.7 Adjust auto shutoff duration


To conserve battery life, the unit automatically turns off after 6 minutes. To customize the auto shutoff setting, press the  (menu) button.


Use the   (arrow) keys to select the Power time menu:



Press the  (menu) button again to enter the value select screen.

Use the   (arrow) keys to select the auto shutoff setting.  
Values: 1–59 minutes or ‘constant on’ (max.)

To return to the main menu screen, press the  (menu) button.

To exit and return to measuring flow, press the  (menu) button again.