

9.0 Troubleshooting

Problem	Possible Cause	Suggested Solution
Sensitivity decreased	Probe clogged Probe line punctured Weak battery	Clean the probe tip to remove any debris Visually inspect probe line for holes* Recharge or return to Restek for battery replacement*
Response decreased	Detector not zeroed	Re-zero detector
LED bar graph stays lit during operation	Detector re-zeroed before unit was purged out Reference gas inlet covered by hand or other object	Allow adequate time for detector to purge, then re-zero Remove obstruction

**Contact Restek or your Restek representative for return instructions for servicing a damaged unit. Additional charges may apply if the warranty has expired or the unit is damaged due to misuse.*