



Figure 2.0

Figure 2.0 shows the complete installation of cables to the GC and for RS 232 remote control. **Be certain all cable connections are made before the power to the Cobra L/S is turned on.**

2.6 Sample Tray Installation

Unwrap the sample tray and place it onto the sample tray brackets. The tray will only mount in one direction as there are alignment pegs on the holder brackets and corresponding alignment slots in the tray. Once installed, place an empty vial with cap and septa in vial position #1 and the waste / solvent locations to be used. Be certain the waste/solvent tray is located on the correct side of the sample tray for your GC. Note, the waste / solvent tray may be located on either side of the sample tray, see Figures 2.6 & 2.6 a.

