



### Cobra L/S Keyboard Overlay

#### Press Hold/Stop Key - One Time.

During any operation of the Auto Sampler the **“Hold/Stop” Key** may be pressed. Hold the key down until the “End Pause” message appears. For an operation pause, the screen will display the prompt **“End Pause”**. The Auto Sampler will continue until the current injection cycle is completed.

At this point practically any aspect of the Auto Sampler may be updated. Methods may be edited, a Priority Manual Sample may be performed or any of the System parameters may be changed. To resume the AutoRun from the point it was paused, press the **“Enter” Key**.

#### Press Hold/Stop Key - Two Times.

Pressing the **“Hold/Stop” Key** the second time will immediately stop the motors. A prompt will appear on the screen to allow the syringe contents to be dumped to the waste vial before returning the syringe to its home position and reset. The method or AutoRun cannot be resumed from the position it was stopped without editing the method and changing the “First Vial Position”. If the syringe contains any solution at this point, press the **“Enter” Key** to allow it to be dumped into the waste vial before returning the syringe to its home position.

**CAUTION:** Syringe contents will be emptied where the syringe is currently positioned if the **“Hold/Stop” key** is pressed instead of the **“Enter” Key**. This could cause hazardous and/or corrosive chemicals to be dispensed into the atmosphere and/or onto the Auto Sampler and Analyzer.