

12	1	Vertical Acceleration	0-255
13	1	Tray Size	0-3 (only 0 currently supported)

5. Cobra will respond with ok.

10.5 System Errors

Undefined commands will get a response "Undefined command"

All parameters must be in their allowed range, Cobra will not edit incoming values.

Cobra will wait for an enter key press (ASCII 13) to continue after an error.

10.6 Error codes

- # 1 Undefined command
- # 2 Invalid value
- # 3 Not yet supported
- # 4 Current run must be aborted first

11.0 Cobra L/S Installation of Cables and Mounting Hardware

The following section details the Cobra L/S mounting bracket hardware and I/O cable installation. Refer to the drawings for detail.

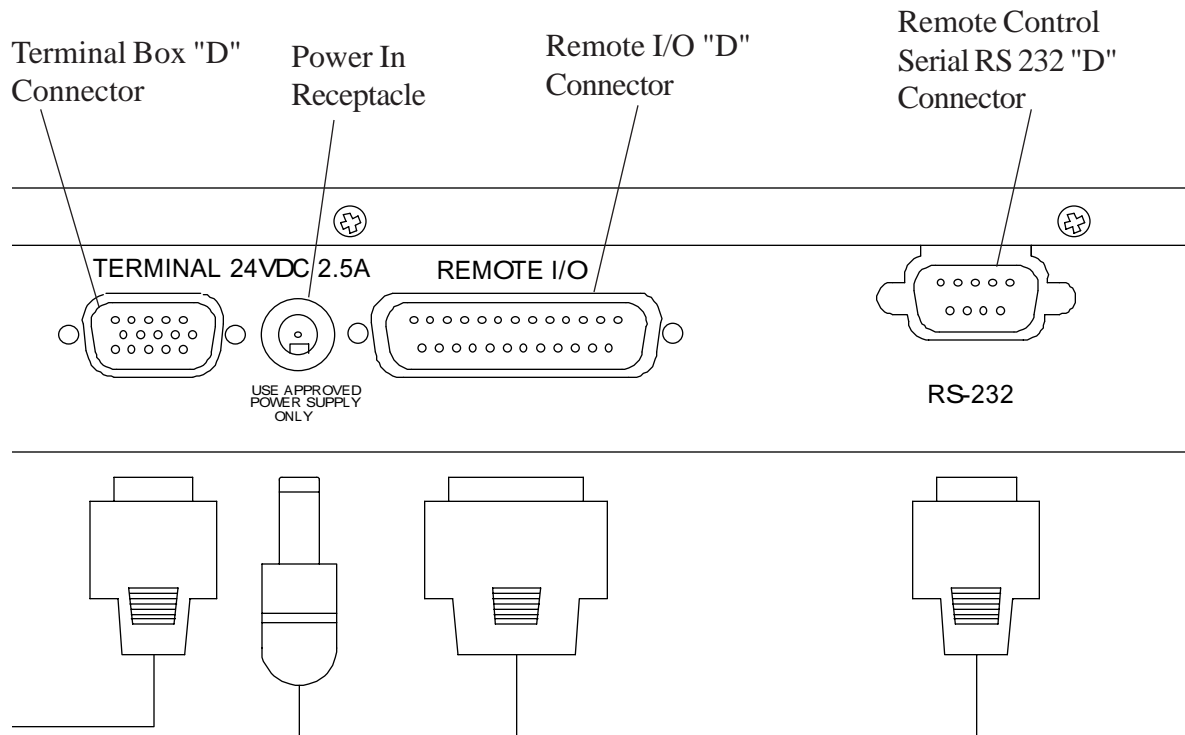


Figure 11.0
Cobra L/S Rear Panel