

Basic Operation of the Manual Remote

The manual remote provides simple valve positioning capabilities, but in the configuration mode, it can be used to perform extensive actuator setup functions. For information on using the manual remote to configure the actuator, refer to "Using the Manual Remote to Configure the Actuator" beginning on the next page.

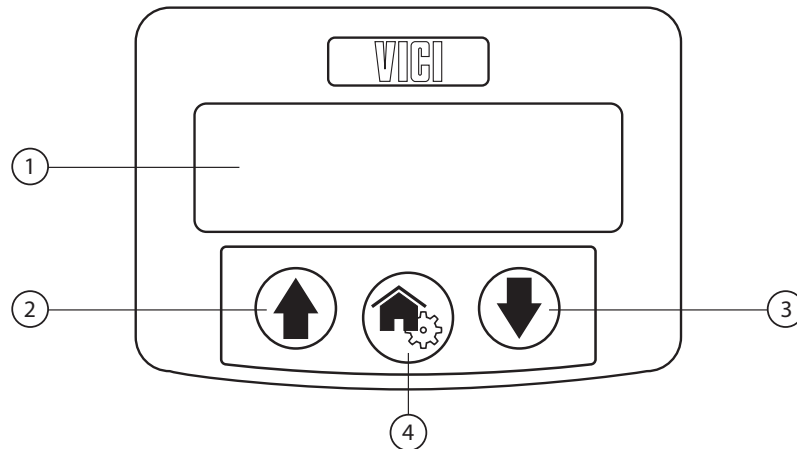


Figure 3: Universal actuator manual remote

1. Display

2. **UP** arrow button

In *two position* configuration, pressing the UP arrow button moves the actuator to Position B. If it is already in Position B, nothing happens.

In *multiposition* configuration, pressing the UP arrow button advances the actuator one position; i.e., from 1 to 2, 4 to 5, etc.

3. **DOWN** arrow button

In *two position* configuration, pressing the DOWN arrow button moves the actuator to Position A. If it is already in Position A, nothing happens.

In *multiposition* configuration, pressing the DOWN arrow button reverses the actuator one position; i.e., from 2 to 1, 5 to 4, etc.

4. **HOME/SETUP** button

In *two position* configuration, pressing the HOME button moves the actuator to Position A. If it is already in Position A, nothing happens.

In *multiposition* configuration, pressing the HOME button sends the actuator to Position 1.

In either configuration, pressing and holding the HOME button for 5 seconds sends the manual remote into the configuration mode. Refer to "Using the Manual Remote to Configure the Actuator" beginning on the next page.