

3. In the Results window locate the **Copy results log** button at the bottom of the window and click on it with the mouse cursor (don't confuse the Copy button with the Copy results log button). Open up Microsoft Excel (or if Excel is not loaded Microsoft Word or Power-Point) and select **Edit** from the menu bar and then **Paste** to copy the results log to Excel.
4. Go back into PeakSimple and close the Results window by selecting the **Close** button. Right click using the mouse cursor on the chromatogram and select **Postrun** from the drop down menu to open the Post-run actions window. From the window locate the Add to results log checkbox and add a check to the box. By selecting the Add to results log checkbox all results from data analysis will automatically be added to the results log after the run is done. Click on **OK** to exit the window. In this way a summary of many analyses can be automatically created and then exported from PeakSimple.



This concludes the PeakSimple 2000 Advanced Tutorial

Further documentation can be obtained by going to:
www.srigc.com online

If you have questions or would like to place an order call:
 (310) 214-5092