

Petrochemical

2011 Applications

RESTEK

2011 • International Distributor Conference



Who are our Customers?

Petroleum



ExxonMobil

Syn crude

Schönbberger

ifp

GAZPROM

Chemicals



WACKER

EASTMAN

elf atochem
ATO

Innophos

RESTEK

2011 • International Distributor Conference



Products analyzed in Segment

Petroleum



Oil, Natural gas, gasoline, kerosene, diesel, Naphtha

Hydrocarbons, Sulfur,
Biomarkers, Aromatics, Alcohols

Bio-Fuels, Biodiesel, Bioethanol

glycerides, FAME, alcohols

Chemicals



Intermediates and final product

C2-C5 Alkenes, Aromatics, Amines,
Glycols, Acids, Solvents, Halogenated
compounds, Alcohols, Aldehydes

RESTEK

2011 • International Distributor Conference



RESTEK

HROM or **ECH** for **ANALYSIS**
ECH for **ANALYSIS**
Website: www.restek.com Email: info@restek.com Tel: 800 676 2016 ... • AUSTRALIA

Australian Distributors
Instruments & Supplies
www.restek.com.au

1213

Outline


- MXT-1HT Sim Dist
- MXT-2887
- Rtx-DHA
- D3606 packed column set
- Rt-XL Sulfur packed column
 - And Siltek deactivation for sulfurs
- PLOT columns
- Appendix
 - Products for biodiesel and chemical analysis

RESTEK

2011 • International Distributor Conference



Refinery Process Map

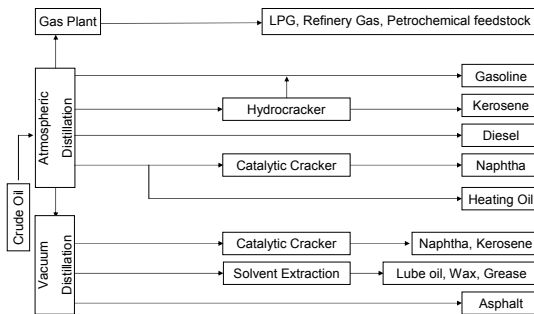

RESTEK

2011 • International Distributor Conference



What does every refinery do?

- Simulated Distillation
- Detailed Hydrocarbon Analysis
- Benzene /Toluene
- Refinery gases
- Sulfur analysis
- Oxygenates

What Solutions
does Restek offer?

RESTEK

2011 • International Distributor Conference




RESTEK

HROM or Zymolene ANALYST 20M
ECH Fresh Air purge for mg dm
Website: www.restek.com Email: info@restek.com Tel: 800 6762 2016 ... • AUSTRALIA

Australian Distributors
Supplies & Instruments
www.restek.com.au

1213

Simulated Distillation

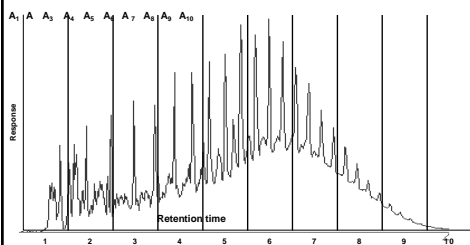

- The chromatogram of the sample is divided into time slices
- The software determines the area count for each slice

RESTEK

2011 • International Distributor Conference



Challenges

- Columns Break (fused silica)
 - Need to recalibrate often (time)
 - Replace columns (price per analysis, time)
 - Reduced accuracy (reliability challenge)
- Columns loose stationary phase by bleed
 - Do not meet resolution requirements

RESTEK

2011 • International Distributor Conference



Needs for Accurate Simulated Distillation Analysis

- High temperature stable columns
- Low Bleed phases

RESTEK

2011 • International Distributor Conference



Sim Dist with fused silica: Big Problem area..

- FS Standard fused silica column polyimide coated, type Inferno..



Polyimide outside coating seriously damaged

This will happen with ALL fused silica columns

RESTEK

2011 • International Distributor Conference



Optimize strength/robustness of capillary Application of MXT Metal capillary tubing

RESTEK

2011 • International Distributor Conference



MXT Tubing is Extremely Strong



- Virtually unbreakable
 - Stainless steel with Siltek coated surface
- Can be bent / coiled to small radius
- Siltek surface stabilizes the stationary phase
- Sim Dist columns wound on standard 11-pin cage
- MXT columns can be cut with standard scoring wafer
 - Special tools also available

MXT columns provide reliability:
No column breakage possible!

RESTEK

2011 • International Distributor Conference



New, easy accessible column cage



MXT is nicely wound on the 11-pin cage:
Easy to install, handle and unwind..



RESTEK

2011 • International Distributor Conference



Simulated Distillation of Crude oil

Analysis: Simulated Distillation

C5-C110 D6352 High Temp

C5-C44 D2887


RESTEK

2011 • International Distributor Conference



ASTM Spec for High Temperature Sim Dist



- In order to be approved for ASTM D86 and 6352, 7169 and 7500:
 - Resolution between C50 and C52 > 2.0, running under ASTM conditions

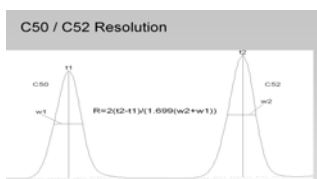

RESTEK

2011 • International Distributor Conference



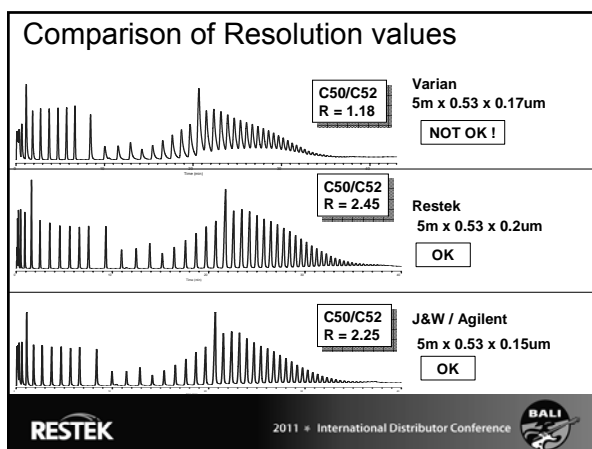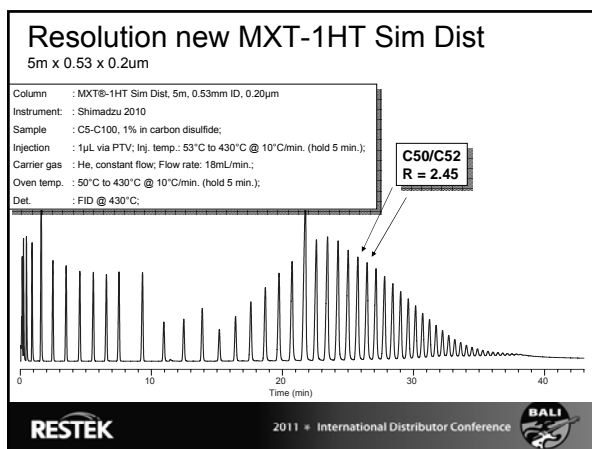

Resolution values

The higher resolution value indicates the Restek MXT-1HT Sim Dist:

- Has a smoother coated surface of the PDMS phase over the length of the column
- The higher resolution value allows for a longer column lifetime

2011 • International Distributor Conference

Column Stability / Lifetime


- Column lifetime in Sim Dist applications are also in direct proportional with the level of the column bleed
- The LOWER the bleed, The LONGER the life time

RESTEK

2011 • International Distributor Conference



Comparison of Bleed values

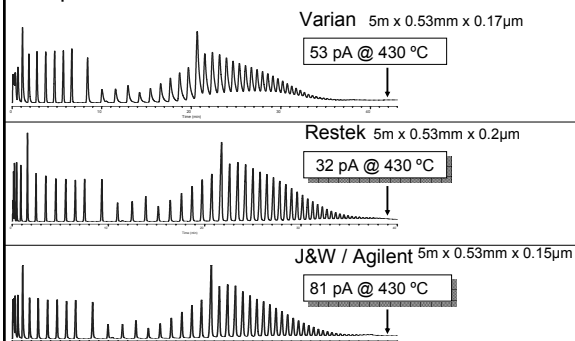

RESTEK

2011 • International Distributor Conference



Performance-ASTM D7169

- Customer feedback on Beta testing, done by Joaquin Lubkowitz, Separation systems.
- Authority in petro – analysis field and ASTM

S Separation Systems, Inc.
Experts In Gas Chromatography

RESTEK



2011 • International Distributor Conference



RESTEK

HROM or Zymol Flow ANALYSIS
ECH Fresh Air purging system
Website: www.separation-systems.com.au Email: info@separation-systems.com.au Tel: 61 07 552 2016 ... • AUSTRALIA

Australian Distributors
Separation Systems, Inc.
1213

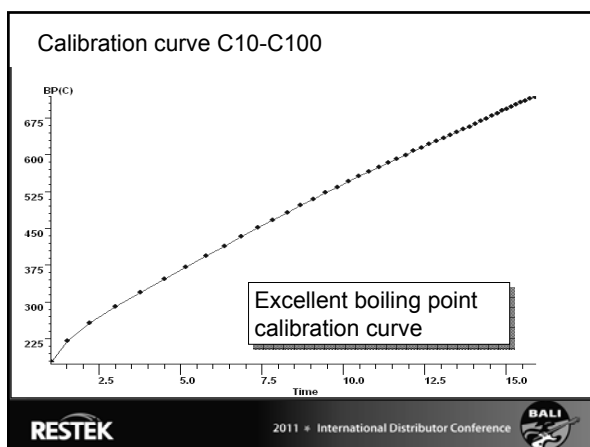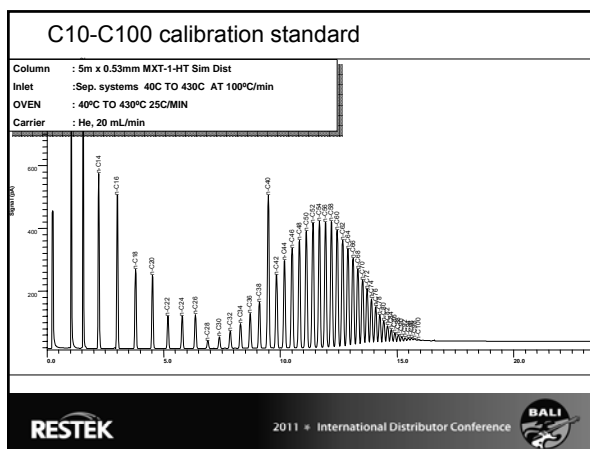

Column Cutting

For installation in a PTV or on-column injector, the column inlet must be well-cut..

RESTEK 2011 • International Distributor Conference BALI

MXT Column Cutting

- A special tube “scorer” has been developed..



Makes a perfect “cut” required for
“on-Column injection”..

For 0.25 - 0.53mm MXT

tubing scorer : 20523
replacement wheel: 20522



RESTEK

2011 • International Distributor Conference



Carrier gas considerations Avoid introduction of oxidizers

RESTEK

2011 • International Distributor Conference



Impact of Oxygen/Moisture

- Increase of bleed
- Retention times drop fast
- Peaks will broaden
- Peaks will tail

RESTEK

2011 • International Distributor Conference



RESTEK

HROM or Zymol Flow ANALYSIS
ECH Fresh Air Supply
Website: www.restek.com.au Email: info@restek.com.au Tel: 02 6762 2016 ... AUSTRALIA


Australian Distributors
Suppliers of Gas Chromatography
www.restek.com.au

1213


Carrier Gas purification

- Degradation reaction of stationary phase + water/oxygen develops exponentially with higher temperature..

Need: Pure gases

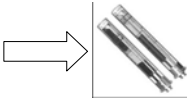

Need: Elimination of leaks

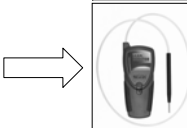

Also use graphite ferrules for temperatures above 360°C



RESTEK

2011 • International Distributor Conference



Simulated Distillation of Gasoline Range Hydrocarbons

Analysis: Simulated Distillation

C5-C110 D6352 High Temp

C5-C44 D2887


Also Called “gasoline-range” Sim Dist

RESTEK

2011 • International Distributor Conference



D2887 Used for:

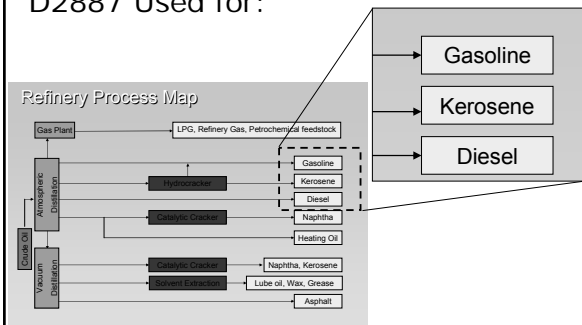

RESTEK

2011 • International Distributor Conference



Gasoline Range Sim Dist

- C5-C44 D2887
- Column: 10m x 0.53mm x 2.65µm Rtx-2887 (Fused silica or MXT)

Uses a THICK film because the sample is injected on the column with no dilution;


Need LOADABILITY !!

RESTEK

2011 • International Distributor Conference



D-2887 Sim Dist

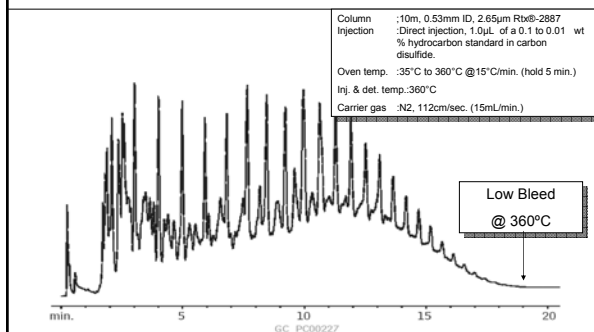

RESTEK

2011 • International Distributor Conference



Restek New Generation MXT Sim Dist Columns

- Features: Low bleeding, Temperature Stable, Better Efficiency
- Benefit: Longer Column Lifetime, Less Replacement, less calibrations, Saving \$\$\$\$\$

RESTEK

2011 • International Distributor Conference



RESTEK

Product information for MXT Sim Dist

- Sim Dist sales companion + presentation
- Restek Refined brochure PCFL1195A: <http://www.restek.com/restek/images/external/PCFL1195A.pdf>
- Simdist brochure PCFL1201A: <http://www.restek.com/restek/images/external/PCFL1201A.pdf>
- Applications see: <http://www.restek.com/restek/class/5004088.asp>; and http://www.restek.com/info_chromatograms.asp
- Technical Literature: http://www.restek.com/aol_lit_main.asp
- Article: Tighten Control of Distillation Processes with the New MXT®-1HT Sim Dist Column : http://www.restek.com/aol_petro_A022.asp



RESTEK

2011 • International Distributor Conference



What does every refinery do?

- Simulated Distillation
- Detailed Hydrocarbon Analysis
- Benzene /Toluene
- Refinery gases
- Sulfur analysis
- Oxygenates

What Solutions
does Restek offer?

RESTEK

2011 • International Distributor Conference



What products come from most refineries?

- Gasoline
- Diesel
- Jet Fuel
- Refinery Gas

RESTEK

2011 • International Distributor Conference



RESTEK



Analysis: Gasoline

Detailed Hydrocarbon Analysis (DHA) D6730

- Column: 100m x 0.25mm x 0.50um Rtx-DHA 100
- Tuning column: 5m x 0.25mm x 1.0um Rtx-5
- Tuning column is for adjusting selectivity for aromatics

RESTEK

2011 • International Distributor Conference



Biggest Challenges for Customers Running ASTM D6730

- Have good peak shape for alcohols
- Have sufficient theoretical plates to separate the complex samples
- Coupling of “tuning” capillary
- Analysis time too long, wish for faster methods
- Reproducibility of column-column quality parameters

RESTEK

2011 • International Distributor Conference



Why do the Refiners Run DHA Analysis?

- At the end a certain gasoline must have a certain octane number..
- It is very important for the refiner to monitor the blending process carefully to avoid too much yield.



Different octane numbers are available

- Accurate DHA analysis can save a lot of money ..



RESTEK

2011 • International Distributor Conference



Important Quality Parameter for DHA Analysis

- The column must not only elute hydrocarbons
- Also need to elute **alcohols**




This is a big challenge and that's why
Rtx-DHA 100 columns are preferred

RESTEK

2011 • International Distributor Conference



Elution of Alcohols on the DHA-type Columns ASTM Method D6730

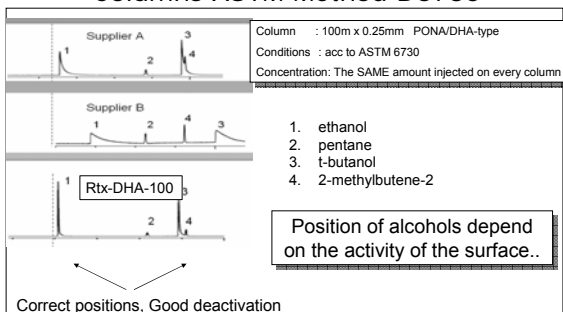

RESTEK

2011 • International Distributor Conference



Impact of a Tailing Alcohol Peak


Supplier 1 type DHA column

Oven: 5 °C → 8.23 min., 22°C/min → 48 °C

24160U

1. ethanol
2. C5
3. 2 methyl butene 2
4. t butanol

Alcohol peaks are retained by column
activity / adsorption;
They elute at wrong position..

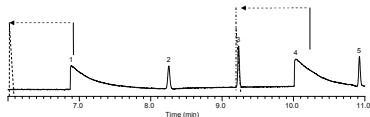

Risk: Misidentification and possible wrong data

RESTEK

2011 • International Distributor Conference



Specifications DHA column: for AstmD6730

Theoretical plates, n, pentane at 35°C;	40000 to 50000
Retention factor, k, pentane at 35°C	0.45 to 0.50
Resolution, R, t-butanol and 2-methylbutene-2 at 35°C	3.25 to 5.25
Peak symmetry, t-butanol at 35°C	>1.0 to <5.0

2011 • International Distributor Conference



Chromatogram showing peaks for ethanol, t-butanol, 2-methyl-butene-2, and pentane. The x-axis is time in minutes (0.00 to 16.00) and the y-axis is signal in pA (-10.00 to 130.00).

Compound	Approximate Retention Time (min)	Approximate Signal (pA)
ethanol	7.5	125
t-butanol	9.5	105
2-methyl-butene-2	9.8	70
pentane	8.8	40

2011 • International Distributor Conference



TEMP PROGRAM: 35°C Isothermal			
INJ TYPE: Split		INJ TEMP: 250°C	
DET TYPE: FID		DET TEMP: 300°C	
LINEAR VELOCITY: 27 cm/sec (Helium)		SPLIT FLOW: 230 mL/min	
Max Cal: 3500			
Retention Time	Name	Area	Height
7.79	ethanol	213288	127.793
9.16	C5	79783	41.430
9.51	1 butanol	226500	108.583
9.80	2 methylbutane 2	162390	77.061
			Width @ 50%
			0.024
			0.028
			0.030
			0.031
Max Cal: 3500			
Partition Ratio (K) C5: 0.48		Resolution (1 butanol / 2 methylbutane 2): 5.63	
Theoretical Plates (C5): 589,096		Asymmetry (1 butanol): 1.38	
Plates per meter (C5): 5,918			
		Column Bleed (pA): 9.9	

2011 • International Distributor Conference




Reproducibility of Rtx-DHA-100 Columns

Following graphs show the typical performance of the chronological production of 250 Rtx-DHA 100 columns

- Efficiency /Resolution
 - Retention
 - Inertness
- } Set by ASTM
- Bleed

RESTEK

2011 • International Distributor Conference

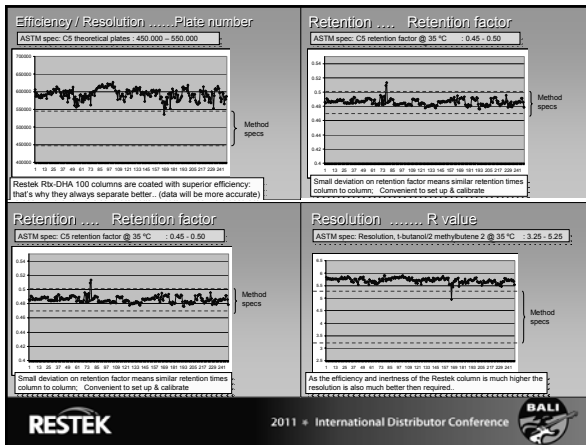

ASTM D 6730 Requires a "tuned" Column...

ASTM wants general purpose methods

Columns from different suppliers will show small differences in selectivity..

By coupling a short piece of Rtx-5, every column can be "tuned" to fulfill ASTM specs on selectivity..




Tuning pre-column: Rtx-5, 2-5 meters
 Length to be determined empirically

RESTEK

2011 • International Distributor Conference



RESTEK

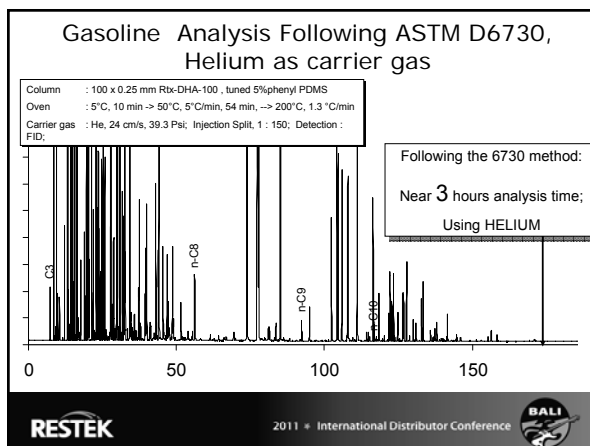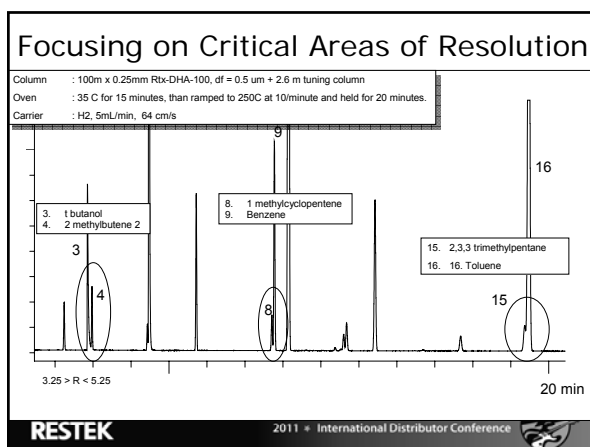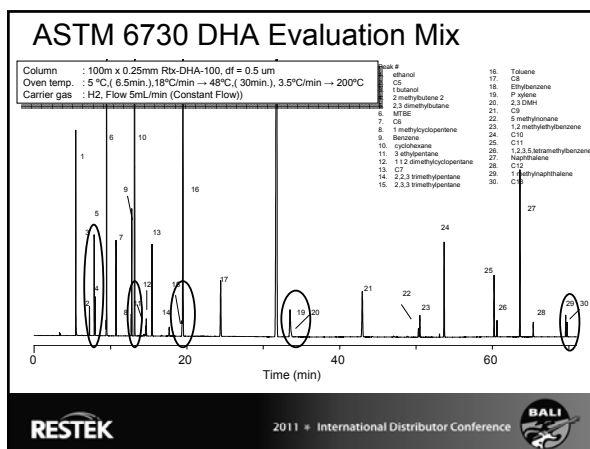

There will be always a wish from customers to SPEED UP the analysis

Using Helium and different temperature program

Using Hydrogen

Using shorter, smaller ID columns



2011 • International Distributor Conference



Using Helium and a different program

Using the same column, same GC

Column : 100m x 0.25mm Rtx-DHA-100, df = 0.5 μ m coupled with
2.6m x 0.25mm Rtx-5DHA tuning column, df = 1.0 μ m

- Carrier gas : He, Flow 2.3 mL/min (Constant Flow), 28 cm/s


- Oven temp. : 5 $^{\circ}$ C \rightarrow 15min., 5 $^{\circ}$ C/min \rightarrow 50 $^{\circ}$ C, 50min.
8 $^{\circ}$ C/min \rightarrow 200 $^{\circ}$ C, 10 min



2011 • International Distributor Conference



Using Helium and a Different Program

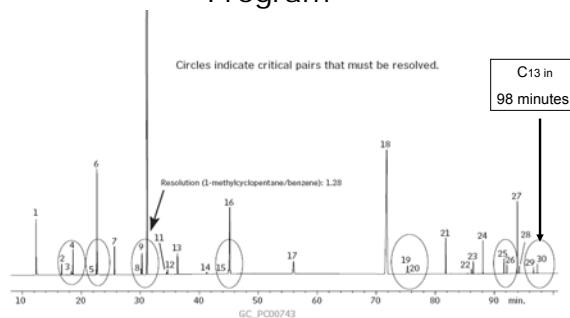

2011 • International Distributor Conference



Speeding up the Application using HYDROGEN Carrier Gas


- Using hydrogen we can operate at near 2 times the linear velocity as used for helium
 - This results in significant shorter run time
- To get the same chromatogram/separation, we need to adjust the oven temperature program

RESTEK

2011 • International Distributor Conference



Fast DHA Analysis using Hydrogen...

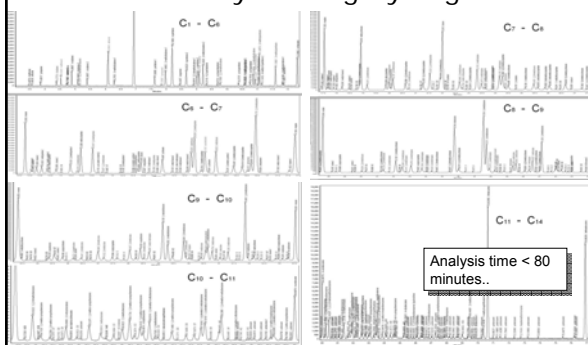

RESTEK

2011 • International Distributor Conference



Fast DHA using Hydrogen

- Same GC
- Same method
- Same Injection/detection techniques
- Same column
- No issues with overloading, peak shifts, tailing, high inlet pressures, discrimination and reduced life time..



Neil Johansen

All retention data and identification is generated by Neil Johansen, developer of the D6730 DHA method.

Available on the Restek website
: http://www.restek.com/promo_gc_pona.asp#004

RESTEK

2011 • International Distributor Conference



RESTEK

HROM or Zymolene ANALYSIS ECH Peak for gas, liquid, and solid samples
Website: www.restek.com Email: info@restek.com Tel: 800 676 2516 ... AUSTRALIA

Australian Distributors
Instruments & Equipment
1213

Hydrogen : Safety Issues

Need to make sure there is no accumulation of H₂ possible in the oven..

Hydrogen is combustible over a concentration range of 4% to 74% by volume;
Diffusion VERY fast;

- Use of Hydrogen generators: limited amount of hydrogen
- Hydrogen Flow restriction to set with GC inlet (electronic flow setting)
- Hydrogen detection systems (sniffer)

RESTEK

2011 • International Distributor Conference



Rtx-DHA Columns for Other ASTM Methods

For ASTM D6733 : Rtx-DHA-50 50m x 0.25mm, df = 0.5 μ m

For ASTM D6729 : Rtx-DHA-100 100m x 0.25mm, df = 0.5 μ m

For ASTM D5501 : Rtx-DHA-150 150m x 0.25mm, df = 1.0 μ m

NOTE:

This column was called earlier Rtx-1 PONA, which is a WRONG name as the application is DHA, not PONA.

Competing products are still called: HP-PONA, DB-PONA, CP-Sil PONA CB, BP-1PONA and Petrocol

RESTEK

2011 • International Distributor Conference



Biggest Challenges for customers running ASTM D6730

Have good peak shape for alcohols.....



Have sufficient theoretical plates



Coupling of "tuning" capillary



Analysis time too long, wish for faster methods ..



Reproducibility of column quality parameters



RESTEK

2011 • International Distributor Conference



RESTEK

HROM or Zymolam HPLC HPLC HPLC
ECH HPLC HPLC HPLC
Website: www.restek.com.au Email: info@restek.com.au Tel: 01 6762 2016 ... • AUSTRALIA

Australian Distributors
Incorporated in Australia
www.restek.com.au

1213

Customer referral

I just wanted to relay to you how pleased I am with the performance of the Restek DHA PONA bonded to the 5 meter RTX-5 pre-column. We've found that the optimal tuning length is 2.05 meters \pm 0.5 meters, which means less time finding the optimal length as compared to the prior Supelco columns.

The peaks look great – far superior to the prior supplier. I am especially pleased with the critical separation of t-butanol and 2-methylbutene-2. The t-butanol peak only tails minimally with the temperature programming, and it keeps us from having to scrap failed columns. This improvement process of switching columns has improved our production rate by nearly half and cut our waste costs as well.

Terry Osenbach
PerkinElmer Health Science
GC Application and Engineered Solutions
Chemist
710 Bridgeport Ave, Shelton, CT 06614

RESTEK

2011 • International Distributor Conference



Product information Rtx DHA columns

- Rtx-DHA sales companion + presentation
- Restek Refined brochure PCFL1195A: <http://www.restek.com/restek/images/external/PCFL1195A.pdf>
Brochure Rtx-DHA 100 PCFL1007A: <http://www.restek.com/restek/images/external/PCFL1007A.pdf>
- Recorded WEBINAR
- Applications see: <http://www.restek.com/restek/class/5004088.asp>; and http://www.restek.com/info_chromatograms.asp
- Technical Literature: http://www.restek.com/aci_lit_main.asp
- Article: **Faster DHA Analyses Using Helium or Hydrogen:** http://www.restek.com/aci_petro_A025.asp
- Article: **How Good is Your PONA Column?:** http://www.restek.com/aci_petro_A006.asp



RESTEK

2011 • International Distributor Conference



What does every refinery do?

- Simulated Distillation
- Detailed Hydrocarbon Analysis
- Benzene / Toluene/ ethanol in gasoline
- Refinery gases
- Sulfur analysis
- Oxygenates

What Solutions
does Restek offer?

RESTEK

2011 • International Distributor Conference




RESTEK



Benzene / Toluene/ Ethanol in Gasoline

ASTM Method D3606

Figure 1. HPLC columns often fail to adequately resolve benzene from ethanol, resulting in poor identification results.

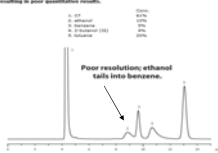

Present solution is problematic:

Big issues with separations and peak shapes

RESTEK

2011 • International Distributor Conference



Modified ASTM D3606 Method using the D3606 Column Set

Uses a SET of packed columns:

Material	Siltek deactivated SS
ID / OD	2mm / 1/8"
Lengths	6' and 16' long
Phases	Rtx-1 and proprietary

- First part of eluting peaks is sent to the second, highly selective column, where the alcohols and aromatics are separated
- Higher boiling compounds are back-flushed
- Unique column set, only from Restek


RESTEK

2011 • International Distributor Conference



New D3606 Method

Columns : Rt™-D3606 Column Set (cat.# 83606)
 column 1 : Rt™-D3606-1: 6'x 1/8"OD (1.8m x 2mm ID), nonpolar RtxB-1
 column 2 : Rt™-D3606-2: 16'x 1/8"OD (4.9m x 2mm ID), proprietary packing
 Injection : 0.05µL, direct injection; Inj. temp. :200°C
 Car. Gas : He, constant flow; Flow rate:20mL/min.
 Oven : 135°C
 Det. : FID, temp: 250°C

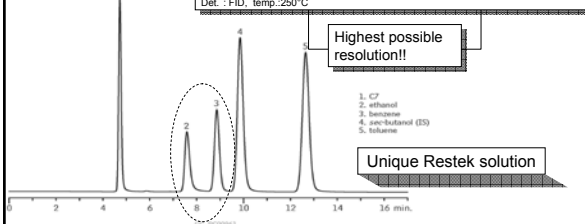

RESTEK

2011 • International Distributor Conference




RESTEK

Australian Distributors
 Repco Pty Ltd
 www.repcopty.com.au

1213

Repeatability of 4 Injections of Walmart Gasoline

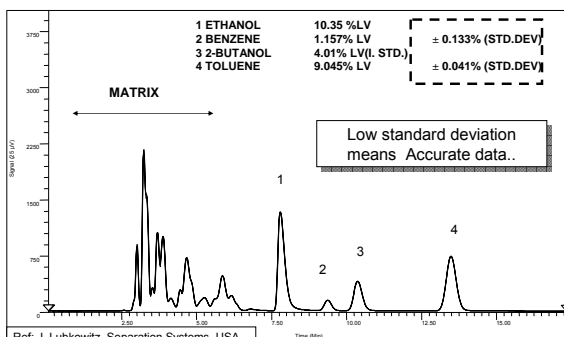

RESTEK

2011 • International Distributor Conference



Product Information for D3606

- Restek Refined brochure: [PCFL1195A: http://www.restek.com/restek/images/external/PCFL1195A.pdf](http://www.restek.com/restek/images/external/PCFL1195A.pdf)
- Applications see : <http://www.restek.com/restek/class/5004088.asp>
- Technical Literature: http://www.restek.com/aoi_lit_main.asp
- Article: **New D3606 Column Set Outperforms TCEP Columns for Benzene Analysis:**
http://www.restek.com/aoi_petro_A020.asp



RESTEK

2011 • International Distributor Conference



What does every refinery do?

- Simulated Distillation
- Detailed Hydrocarbon Analysis
- Benzene / Toluene/ ethanol in gasoline
- Refinery gases
- Sulfur analysis
- Oxygenates

What Solutions
does Restek offer?

RESTEK

2011 • International Distributor Conference



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
RESTEK




PLOT Layer is often deposited as non homogeneous film

FLOW Restriction

Ideal "open Path: Flow-restriction factor = 1.0



PLOT column with Flow-Restrictions: Flow restriction factor "F" = 0.2-0.8

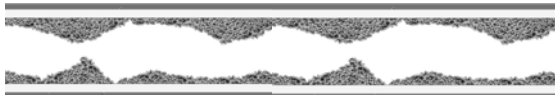

RESTEK

2011 • International Distributor Conference



This was one of the reasons why we could replace Varian and Agilent PLOT columns for:

- Siemens
- Agilent micro GC
- C2V (Thermo)
- Schlumberger
- Many companies that supply Analyzer solutions (PAC, Interscience,...)

RESTEK

2011 • International Distributor Conference




Transportation



Transportation of PLOT columns

During transport columns are subjected to uncontrolled:

- Vibrations
- Pressure changes
- Temperature changes
- Shocks

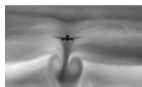

Which all can destabilize the adsorption layer..

RESTEK

2011 • International Distributor Conference



RESTEK

HROM or Zymol Flow 40/60 100/20 150/30
ECH Flow 40/60 100/20 150/30
Website: www.restek.com.au Email: info@restek.com.au Tel: 61 6762 2016 ... @ AUSTRALIA

Australian Distributors
Suppliers of Gas Chromatography
www.restek.com.au

12/13

Dow Chemical, J . Luong

One of the concerns involves the use of PLOT column technology is adsorbent particle shedding. The shedding of adsorbent particles has the potential of plugging the flow channel of the capillary flow Deans switch device. This in turn can cause retention time shift.

Prior to connecting the columns to the capillary flow Deans switch device, we highly recommend the columns be purged with carrier gas at a flow of 50 mL/min for 15 min to remove any particles that might have been dislodged during shipment of the columns from the manufacturers.



2011 • International Distributor Conference



Refinery Gas Analysis

- C1-C5
 - Rt AluminaBOND KCl or Na2SO4
- Permanent gases, C1-C2 and CO2:
 - Rt Msieve 5A / Rt-Q BOND
 - ShinCarbon



2011 • International Distributor Conference



C1-C5 Hydrocarbons Rt-Alumina BOND

KCl : Non polar Al2O3 surface




Na2SO4 : Polar Al2O3 surface




Deactivation is necessary to reduce the activity of the alumina..



2011 • International Distributor Conference




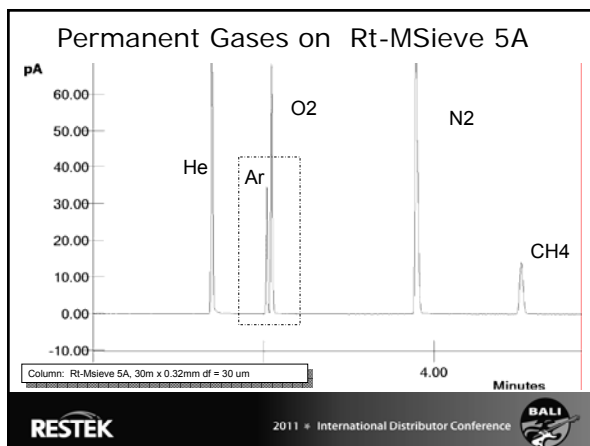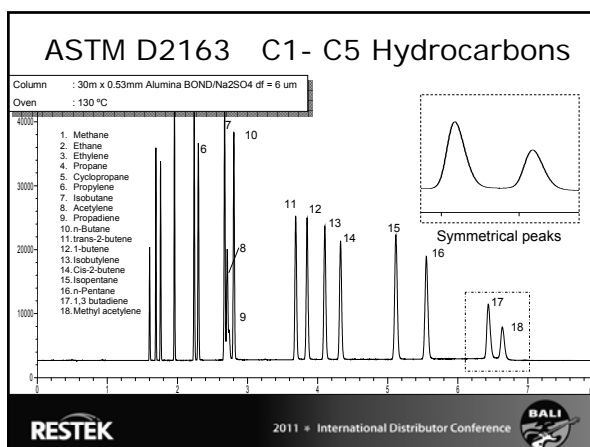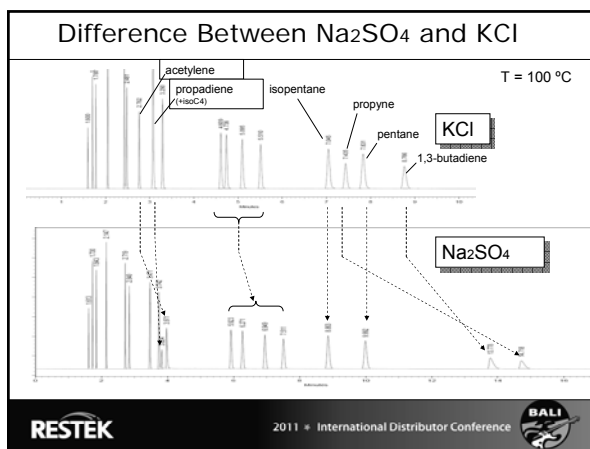

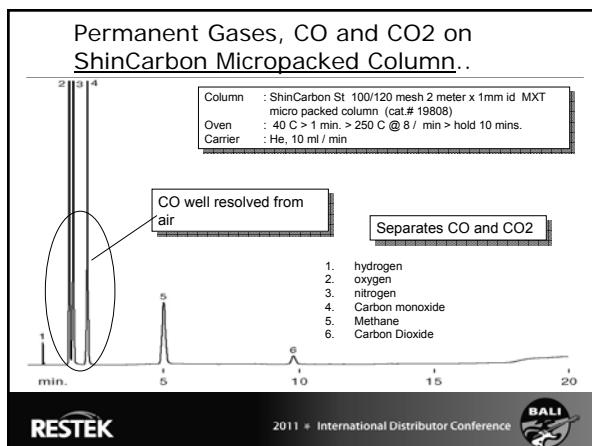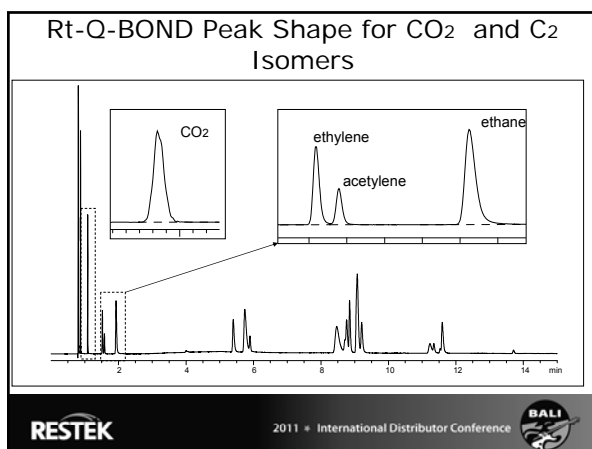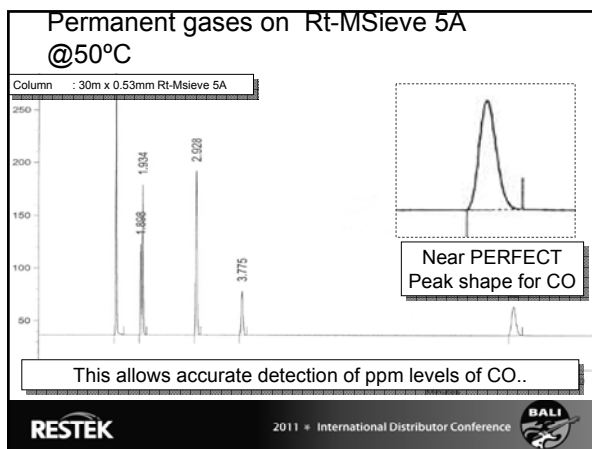

Product Information for PLOT

- Restek Refined brochure PCFL1195A: <http://www.restek.com/restek/images/external/PCFL1195A.pdf>
- Next generation PLOT columns PCFL1163A: <http://www.restek.com/restek/images/external/PCFL1163A.pdf>
- Applications see: <http://www.restek.com/restek/class/5004088.asp>
- Technical Literature: http://www.restek.com/aol_lit_main.asp
- Article: New PLOT Column Technology Stabilizes Flow and Improves Analysis of Light Hydrocarbons, Solvents, and Permanent Gases: http://www.restek.com/aol_petro_A023.asp
- Article: Advances in Porous Layer Open Tubular Columns: http://www.icpubs.com/Media/PublishingTitles/dec/eeuw%20layout_0.pdf

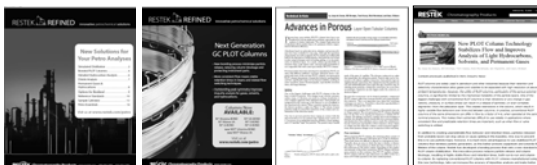

RESTEK

2011 • International Distributor Conference



Additional UNIQUE Restek Development

Making PLOT and LIQUID PHASE columns available in MXT..

In PROCESS type applications, customers look for:

RELIABILITY..

MXT columns will offer extra security and value..

RESTEK

2011 • International Distributor Conference



MXT-MSieve 5A



RESTEK

2011 • International Distributor Conference






RESTEK




HROM or Zymolam 40/60 50/60 50/80 50/100 50/120 50/150 50/200 50/250 50/300 50/350 50/400 50/450 50/500 50/550 50/600 50/650 50/700 50/750 50/800 50/850 50/900 50/950 50/1000 50/1100 50/1200 50/1300 50/1400 50/1500 50/1600 50/1700 50/1800 50/1900 50/2000 50/2100 50/2200 50/2300 50/2400 50/2500 50/2600 50/2700 50/2800 50/2900 50/3000 50/3100 50/3200 50/3300 50/3400 50/3500 50/3600 50/3700 50/3800 50/3900 50/4000 50/4100 50/4200 50/4300 50/4400 50/4500 50/4600 50/4700 50/4800 50/4900 50/5000 50/5100 50/5200 50/5300 50/5400 50/5500 50/5600 50/5700 50/5800 50/5900 50/6000 50/6100 50/6200 50/6300 50/6400 50/6500 50/6600 50/6700 50/6800 50/6900 50/7000 50/7100 50/7200 50/7300 50/7400 50/7500 50/7600 50/7700 50/7800 50/7900 50/8000 50/8100 50/8200 50/8300 50/8400 50/8500 50/8600 50/8700 50/8800 50/8900 50/9000 50/9100 50/9200 50/9300 50/9400 50/9500 50/9600 50/9700 50/9800 50/9900 50/10000 50/10100 50/10200 50/10300 50/10400 50/10500 50/10600 50/10700 50/10800 50/10900 50/11000 50/11100 50/11200 50/11300 50/11400 50/11500 50/11600 50/11700 50/11800 50/11900 50/12000 50/12100 50/12200 50/12300 50/12400 50/12500 50/12600 50/12700 50/12800 50/12900 50/13000 50/13100 50/13200 50/13300 50/13400 50/13500 50/13600 50/13700 50/13800 50/13900 50/14000 50/14100 50/14200 50/14300 50/14400 50/14500 50/14600 50/14700 50/14800 50/14900 50/15000 50/15100 50/15200 50/15300 50/15400 50/15500 50/15600 50/15700 50/15800 50/15900 50/16000 50/16100 50/16200 50/16300 50/16400 50/16500 50/16600 50/16700 50/16800 50/16900 50/17000 50/17100 50/17200 50/17300 50/17400 50/17500 50/17600 50/17700 50/17800 50/17900 50/18000 50/18100 50/18200 50/18300 50/18400 50/18500 50/18600 50/18700 50/18800 50/18900 50/19000 50/19100 50/19200 50/19300 50/19400 50/19500 50/19600 50/19700 50/19800 50/19900 50/20000 50/20100 50/20200 50/20300 50/20400 50/20500 50/20600 50/20700 50/20800 50/20900 50/21000 50/21100 50/21200 50/21300 50/21400 50/21500 50/21600 50/21700 50/21800 50/21900 50/22000 50/22100 50/22200 50/22300 50/22400 50/22500 50/22600 50/22700 50/22800 50/22900 50/23000 50/23100 50/23200 50/23300 50/23400 50/23500 50/23600 50/23700 50/23800 50/23900 50/24000 50/24100 50/24200 50/24300 50/24400 50/24500 50/24600 50/24700 50/24800 50/24900 50/25000 50/25100 50/25200 50/25300 50/25400 50/25500 50/25600 50/25700 50/25800 50/25900 50/26000 50/26100 50/26200 50/26300 50/26400 50/26500 50/26600 50/26700 50/26800 50/26900 50/27000 50/27100 50/27200 50/27300 50/27400 50/27500 50/27600 50/27700 50/27800 50/27900 50/28000 50/28100 50/28200 50/28300 50/28400 50/28500 50/28600 50/28700 50/28800 50/28900 50/29000 50/29100 50/29200 50/29300 50/29400 50/29500 50/29600 50/29700 50/29800 50/29900 50/30000 50/30100 50/30200 50/30300 50/30400 50/30500 50/30600 50/30700 50/30800 50/30900 50/31000 50/31100 50/31200 50/31300 50/31400 50/31500 50/31600 50/31700 50/31800 50/31900 50/32000 50/32100 50/32200 50/32300 50/32400 50/32500 50/32600 50/32700 50/32800 50/32900 50/33000 50/33100 50/33200 50/33300 50/33400 50/33500 50/33600 50/33700 50/33800 50/33900 50/34000 50/34100 50/34200 50/34300 50/34400 50/34500 50/34600 50/34700 50/34800 50/34900 50/35000 50/35100 50/35200 50/35300 50/35400 50/35500 50/35600 50/35700 50/35800 50/35900 50/36000 50/36100 50/36200 50/36300 50/36400 50/36500 50/36600 50/36700 50/36800 50/36900 50/37000 50/37100 50/37200 50/37300 50/37400 50/37500 50/37600 50/37700 50/37800 50/37900 50/38000 50/38100 50/38200 50/38300 50/38400 50/38500 50/38600 50/38700 50/38800 50/38900 50/39000 50/39100 50/39200 50/39300 50/39400 50/39500 50/39600 50/39700 50/39800 50/39900 50/40000 50/40100 50/40200 50/40300 50/40400 50/40500 50/40600 50/40700 50/40800 50/40900 50/41000 50/41100 50/41200 50/41300 50/41400 50/41500 50/41600 50/41700 50/41800 50/41900 50/42000 50/42100 50/42200 50/42300 50/42400 50/42500 50/42600 50/42700 50/42800 50/42900 50/43000 50/43100 50/43200 50/43300 50/43400 50/43500 50/43600 50/43700 50/43800 50/43900 50/44000 50/44100 50/44200 50/44300 50/44400 50/44500 50/44600 50/44700 50/44800 50/44900 50/45000 50/45100 50/45200 50/45300 50/45400 50/45500 50/45600 50/45700 50/45800 50/45900 50/46000 50/46100 50/46200 50/46300 50/46400 50/46500 50/46600 50/46700 50/46800 50/46900 50/47000 50/47100 50/47200 50/47300 50/47400 50/47500 50/47600 50/47700 50/47800 50/47900 50/48000 50/48100 50/48200 50/48300 50/48400 50/48500 50/48600 50/48700 50/48800 50/48900 50/49000 50/49100 50/49200 50/49300 50/49400 50/49500 50/49600 50/49700 50/49800 50/49900 50/50000 50/50100 50/50200 50/50300 50/50400 50/50500 50/50600 50/50700 50/50800 50/50900 50/51000 50/51100 50/51200 50/51300 50/51400 50/51500 50/51600 50/51700 50/51800 50/51900 50/52000 50/52100 50/52200 50/52300 50/52400 50/52500 50/52600 50/52700 50/52800 50/52900 50/53000 50/53100 50/53200 50/53300 50/53400 50/53500 50/53600 50/53700 50/53800 50/53900 50/54000 50/54100 50/54200 50/54300 50/54400 50/54500 50/54600 50/54700 50/54800 50/54900 50/55000 50/55100 50/55200 50/55300 50/55400 50/55500 50/55600 50/55700 50/55800 50/55900 50/56000 50/56100 50/56200 50/56300 50/56400 50/56500 50/56600 50/56700 50/56800 50/56900 50/57000 50/57100 50/57200 50/57300 50/57400 50/57500 50/57600 50/57700 50/57800 50/57900 50/58000 50/58100 50/58200 50/58300 50/58400 50/58500 50/58600 50/58700 50/58800 50/58900 50/59000 50/59100 50/59200 50/59300 50/59400 50/59500 50/59600 50/59700 50/59800 50/59900 50/60000 50/60100 50/60200 50/60300 50/60400 50/60500 50/60600 50/60700 50/60800 50/60900 50/61000 50/61100 50/61200 50/61300 50/61400 50/61500 50/61600 50/61700 50/61800 50/61900 50/62000 50/62100 50/62200 50/62300 50/62400 50/62500 50/62600 50/62700 50/62800 50/62900 50/63000 50/63100 50/63200 50/63300 50/63400 50/63500 50/63600 50/63700 50/63800 50/63900 50/64000 50/64100 50/64200 50/64300 50/64400 50/64500 50/64600 50/64700 50/64800 50/64900 50/65000 50/65100 50/65200 50/65300 50/65400 50/65500 50/65600 50/65700 50/65800 50/65900 50/66000 50/66100 50/66200 50/66300 50/66400 50/66500 50/66600 50/66700 50/66800 50/66900 50/67000 50/67100 50/67200 50/67300 50/67400 50/67500 50/67600 50/67700 50/67800 50/67900 50/68000 50/68100 50/68200 50/68300 50/68400 50/68500 50/68600 50/68700 50/68800 50/68900 50/69000 50/69100 50/69200 50/69300 50/69400 50/69500 50/69600 50/69700 50/69800 50/69900 50/70000 50/70100 50/70200 50/70300 50/70400 50/70500 50/70600 50/70700 50/70800 50/70900 50/71000 50/71100 50/71200 50/71300 50/71400 50/71500 50/71600 50/71700 50/71800 50/71900 50/72000 50/72100 50/72200 50/72300 50/72400 50/72500 50/72600 50/72700 50/72800 50/72900 50/73000 50/73100 50/73200 50/73300 50/73400 50/73500 50/73600 50/73700 50/73800 50/73900 50/74000 50/74100 50/74200 50/74300 50/74400 50/74500 50/74600 50/74700 50/74800 50/74900 50/75000 50/75100 50/75200 50/75300 50/75400 50/75500 50/75600 50/75700 50/75800 50/75900 50/76000 50/76100 50/76200 50/76300 50/76400 50/76500 50/76600 50/76700 50/76800 50/76900 50/77000 50/77100 50/77200 50/77300 50/77400 50/77500 50/77600 50/77700 50/77800 50/77900 50/78000 50/78100 50/78200 50/78300 50/78400 50/78500 50/78600 50/78700 50/78800 50/78900 50/79000 50/79100 50/79200 50/79300 50/79400 50/79500 50/79600 50/79700 50/79800 50/79900 50/80000 50/80100 50/80200 50/80300 50/80400 50/80500 50/80600 50/80700 50/80800 50/80900 50/81000 50/81100 50/81200 50/81300 50/81400 50/81500 50/81600 50/81700 50/81800 50/81900 50/82000 50/82100 50/82200 50/82300 50/82400 50/82500 50/82600 50/82700 50/82800 50/82900 50/83000 50/83100 50/83200 50/83300 50/83400 50/83500 50/83600 50/83700 50/83800 50/83900 50/84000 50/84100 50/84200 50/84300 50/84400 50/84500 50/84600 50/84700 50/84800 50/84900 50/85000 50/85100 50/85200 50/85300 50/85400 50/85500 50/85600 50/85700 50/85800 50/85900 50/86000 50/86100 50/86200 50/86300 50/86400 50/86500 50/86600 50/86700 50/86800 50/86900 50/87000 50/87100 50/87200 50/87300 50/87400 50/87500 50/87600 50/87700 50/87800 50/87900 50/88000 50/88100 50/88200 50/88300 50/88400 50/88500 50/88600 50/88700 50/88800 50/88900 50/89000 50/89100 50/89200 50/89300 50/89400 50/89500 50/89600 50/89700 50/89800 50/89900 50/90000 50/90100 50/90200 50/90300 50/90400 50/90500 50/90600 50/90700 50/90800 50/90900 50/91000 50/91100 50/91200 50/91300 50/91400 50/91500 50/91600 50/91700 50/91800 50/91900 50/92000 50/92100 50/92200 50/92300 50/92400 50/92500 50/92600 50/92700 50/92800 50/92900 50/93000 50/93100 50/93200 50/93300 50/93400 50/93500 50/93600 50/93700 50/93800 50/93900 50/94000 50/94100 50/94200 50/94300 50/94400 50/94500 50/94600 50/94700 50/94800 50/94900 50/95000 50/95100 50/95200 50/95300 50/95400 50/95500 50/95600 50/95700 50/95800 50/95900 50/96000 50/96100 50/96200 50/96300 50/96400 50/96500 50/96600 50/96700 50/96800 50/96900 50/97000 50/97100 50/97200 50/97300 50/97400 50/97500 50/97600 50/97700 50/97800 50/97900 50/98000 50/98100 50/98200 50/98300 50/98400 50/98500 50/98600 50/98700 50/98800 50/98900 50/99000 50/99100 50/99200 50/99300 50/99400 50/99500 50/99600 50/99700 50/99800 50/99900 50/100000 50/100100 50/100200 50/100300 50/100400 50/100500 50/100600 50/100700 50/100800 50/100900 50/101000 50/101100 50/101200 50/101300 50/101400 50/101500 50/101600 50/101700 50/101800 50/101900 50/102000 50/102100 50/102200 50/102300 50/102400 50/102500 50/102600 50/102700 50/102800 50/102900 50/103000 50/103100 50/103200 50/103300 50/103400 50/103500 50/103600 50/103700 50/103800 50/103900 50/104000 50/104100 50/104200 50/104300 50/104400 50/104500 50/104600 50/104700 50/104800 50/104900 50/105000 50/105100 50/105200 50/105300 50/105400 50/105500 50/105600 50/105700 50/105800 50/105900 50/106000 50/106100 50/106200 50/106300 50/106400 50/106500 50/106600 50/106700 50/106800 50/106900 50/107000 50/107100 50/107200 50/107300 50/107400 50/107500 50/107600 50/107700 50/107800 50/107900 50/108000 50/108100 50/108200 50/108300 50/108400 50/108500 50/108600 50/108700 50/108800 50/108900 50/109000 50/109100 50/109200 50/109300 50/109400 50/109500 50/109600 50/109700 50/109800 50/109900 50/110000 50/110100 50/110200 50/110300 50/110400 50/110500 50/110600 50/110700 50/110800 50/110900 50/111000 50/111100 50/111200 50/111300 50/111400 50/111500 50/111600 50/111700 50/111800 50/111900 50/112000 50/112100 50/112200 50/112300 50/112400 50/112500 50/112600 50/112700 50/112800 50/112900 50/113000 50/113100 50/113200 50/113300 50/113400 50/113500 50/113600 50/113700 50/113800 50/113900 50/114000 50/114100 50/114200 50/114300 50/114400 50/114500 50/114600 50/114700 50/114800 50/114900 50/115000 50/115100 50/115200 50/115300 50/115400 50/115500 50/115600 50/115700 50/115800 50/115900 50/116000 50/116100 50/116200 50/116300 50/116400 50/116500 50/116600 50/116700 50/116800 50/116900 50/117000 50/117100 50/117200 50/117300 50/117400 50/117500 50/117600 50/117700 50/117800 50/117900 50/118000 50/118100 50/118200 50/118300 50/118400 50/118500 50/118600 50/118700 50/118800 50/118900 50/119000 50/119100 50/119200 50/119300 50/119400 50/119500 50/119600 50/119700 50/119800 50/119900 50/120000 50/120100 50/120200 50/120300 50/120400 50/120500 50/120600 50/120700 50/120800 50/120900 50/121000 50/121100 50/121200 50/121300 50/121400 50/121500 50/121600 50/121700 50/12

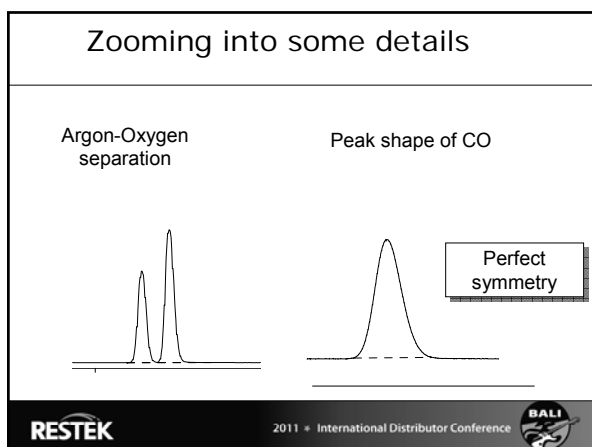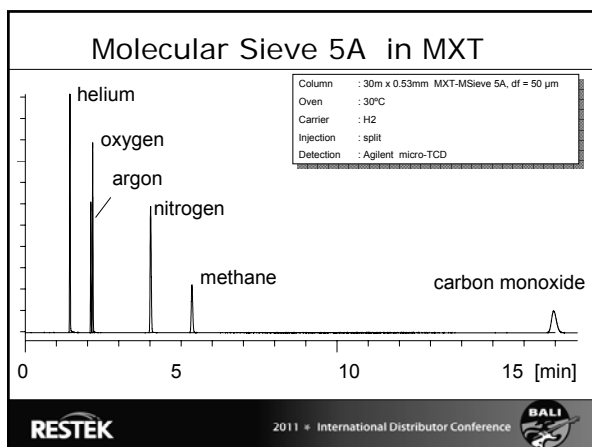

MXT-Alumina BOND (Na₂SO₄ deactivation)

RESTEK 2011 • International Distributor Conference

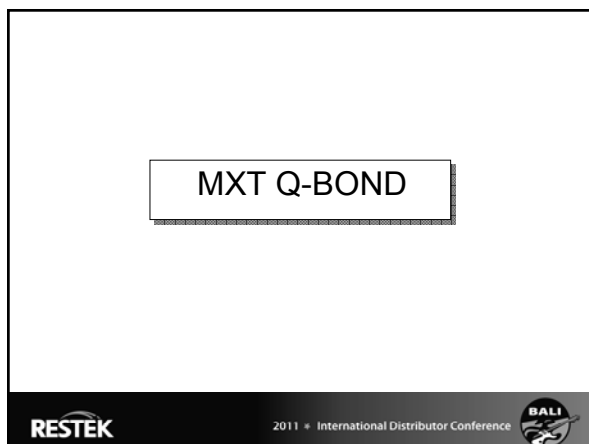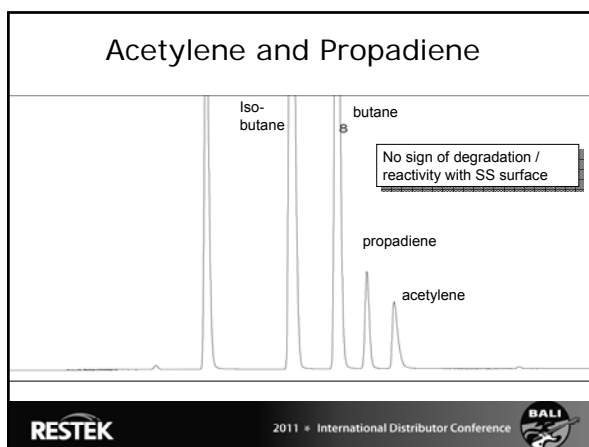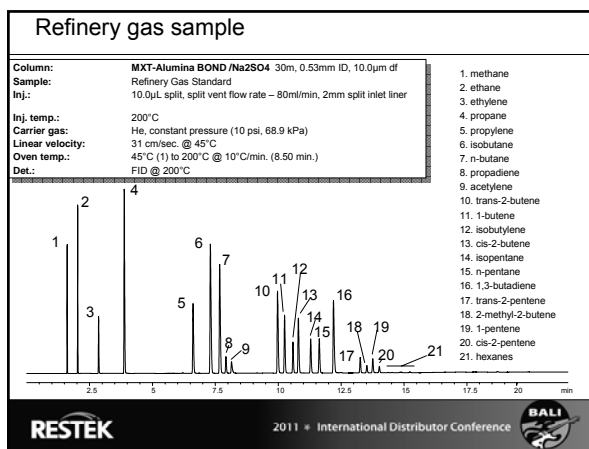

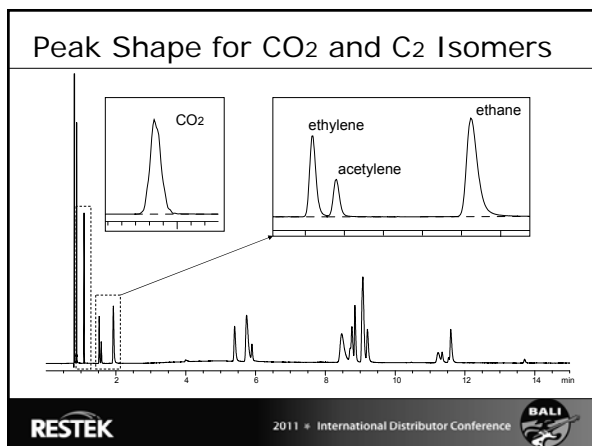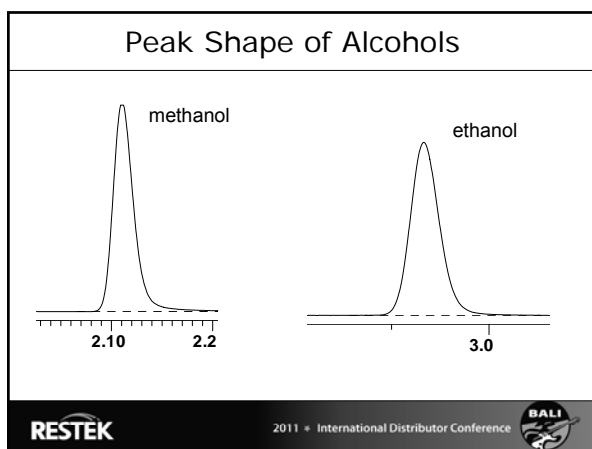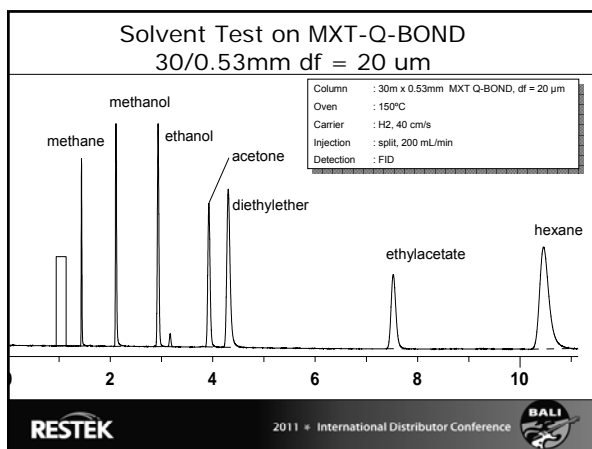

Product information MXT PLOT

- Brochure Restek Refined PCFL1195A: <http://www.restek.com/restek/images/external/PCFL1195A.pdf>
- Brochure Next generation PLOT columns PCFL1163A: <http://www.restek.com/restek/images/external/PCFL1163A.pdf>
- Applications see: <http://www.restek.com/restek/class/5004088.asp> and http://www.restek.com/info_chromatograms.asp
- Technical Literature: http://www.restek.com/aci_lit_main.asp



RESTEK

2011 • International Distributor Conference



What does every refinery do?

- Simulated Distillation
- Detailed Hydrocarbon Analysis
- Benzene / Toluene/ ethanol in gasoline
- Refinery gases
- Sulfur analysis
- Oxygenates

What Solutions
does Restek offer?

RESTEK

2011 • International Distributor Conference



Reference

Restek is internationally known for the Siltek / and SilcoSteel deactivations;


Siltek treated parts are widely used and acknowledged in the petrochemical Industry for sulfur compound stability...

RESTEK

2011 • International Distributor Conference



All parts in the GC system that come in contact with sulfur compounds can be deactivated by the Siltek technology...

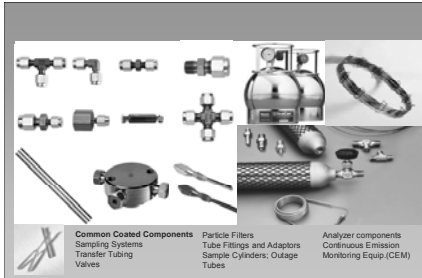


RESTEK

2011 • International Distributor Conference



Sulfur Gases at 19 ppb

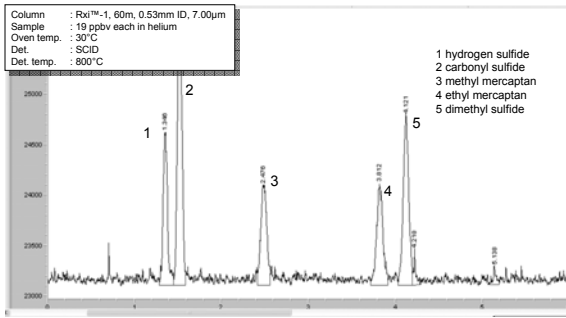

RESTEK

2011 • International Distributor Conference



Most Important Issue for Trace Sulfur Analysis

System Inertness

Use a complete Inert system

- Deactivated transfer tubing, Inert Valves / coupling devices, Inert columns, Inert detection liners, Pressure regulators




Sulfur compounds will disappear especially at low levels..

Restek Developed Siltek deactivated tubing..

RESTEK

2011 • International Distributor Conference



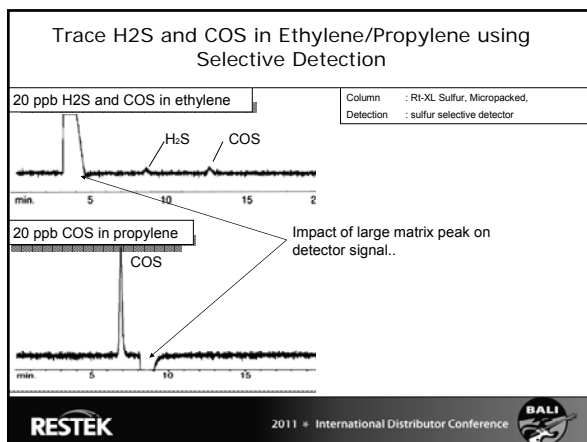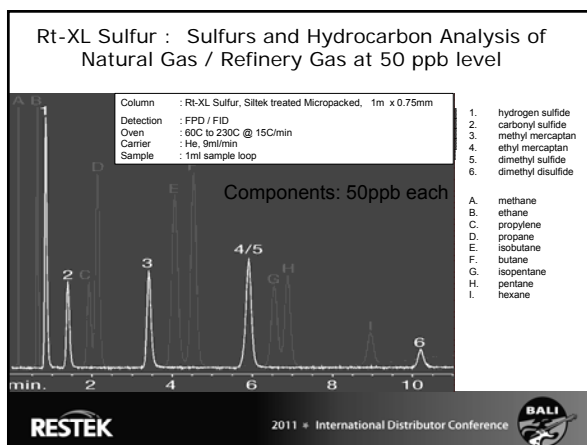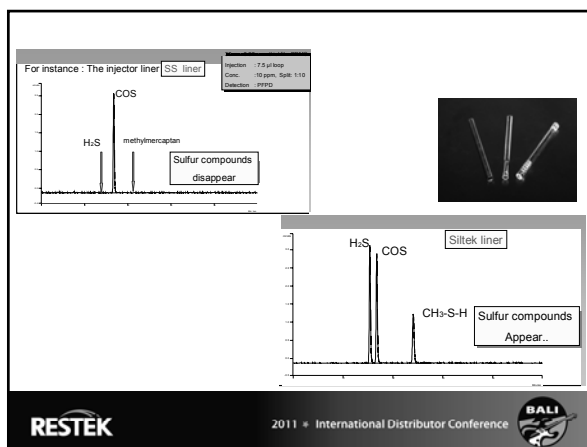

Product information Sulfur

- Restek Refined brochure : PCFL1195A; <http://www.restek.com/restek/images/external/PCFL1195A.pdf>
- Applications : [gc_pc00202.pdf](#); [gc_pc00202.pdf](#); [gc_pc00202.pdf](#); [gc_pc00199.pdf](#); [gc_pc00199.pdf](#)
- Powerpoint handout : http://www.restek.com/restek/images/external/PASS6_SF_TraceLevelSulfurs_Restek.pdf
- Applications see : <http://www.restek.com/restek/class/504088.asp>
- Technical Literature : http://www.restek.com/aol_111_main.asp
- Article: Stable Sulfur & Mercury Sampling in Refineries Using Siltek® and Sulfinet® Surface Treated Components: http://www.restek.com/aol_petro_A016.asp



RESTEK

2011 • International Distributor Conference



What does every refinery do?

- Simulated Distillation
- Detailed Hydrocarbon Analysis
- Benzene / Toluene/ ethanol in gasoline
- Refinery gases
- Sulfur analysis
- Oxygenates

What Solutions
does Restek offer?

RESTEK

2011 • International Distributor Conference



Oxygenates in Gasoline Analyses: ASTM D4815

- Oxygenates in Finished Gasoline (Octane boosters) D4815
- Determination of MTBE, ETBE, TAME, DIPE, tertiary-Amyl Alcohol and C1 to C4 Alcohols in Gasoline by Gas Chromatography
- TCEP column plus Rtx-1 column or Rtx-DHA column

RESTEK

2011 • International Distributor Conference




RESTEK

HROM or Zymolene 95/5 95/5 95/5 95/5
ECH 100% 100% 100% 100%
Website: www.restek.com Email: info@restek.com Tel: 800 6762 2016 ... in AUSTRALIA

Australian Distributors
Instruments & Electronics
www.restek.com.au

1213

Alcohols in Gasoline by ASTM D4815

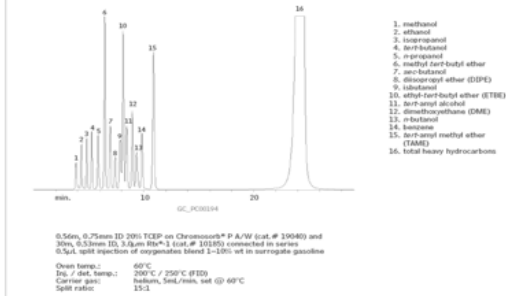

*Silcosteel® treatment is a proprietary surface treatment for passivating steel and stainless steel, U.S. Patent 6,513,760.


RESTEK

2011 • International Distributor Conference



Product Information for Oxygenates

- Applications : gc_pc00194.pdf
- Application article : Oxygenates in gasoline: 59587A

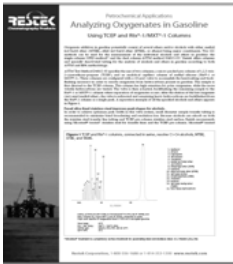

RESTEK

2011 • International Distributor Conference



RESTEK

Website: www.restek.com Email: info@restek.com Tel: 800 678 2636 ... @ RESTEK